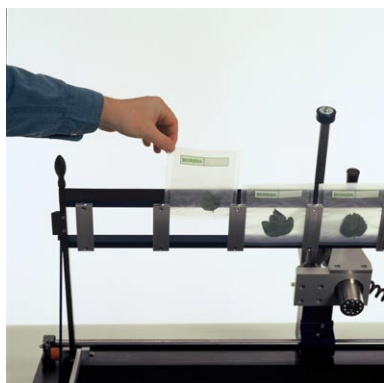
Product Information

HOMEX 6 Instructions

Filling of the rack

Insert 1 to 5 extraction bags in the open rack.

Closing the rack

Close the rack by turning the two clamping-screws (1) to fix the bags in the homogenization position.

Homogenization

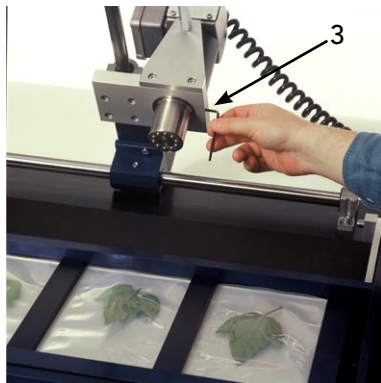
Switch the motor on and homogenize the sample by gliding the „ball-bearing head“ over the bags, using the handle.

Changing the rack


To change the rack, push the knobs (2) at the two fixing-screws and pull the screws out. Insert the rack and put the screws back in position.

Removing the ball-bearing head

To change or clean the ball-bearing head, open first the screw on the side (3) and remove the cover-case. To remove the ball-bearing head entirely, the head fixation screw (4) has to be removed as well.

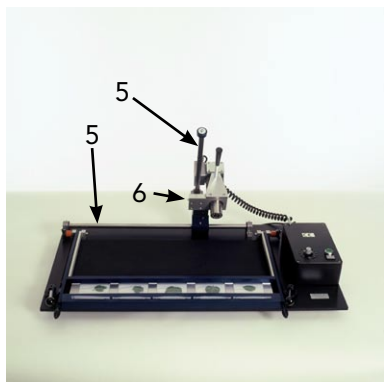Maintenance and storage

Remove spills of extraction buffer immediately first with a soft cloth wetted with water, then wipe dry - in particular when buffer entered the ball-bearing head.

Occasionally grease the two gliding axes (5) and the nipple (6) at the homogenizing unit. Check the holding-screws of the motor and the rack and tighten them if necessary. Replace the aluminium bars holding the bags when they look worn. They are made of „soft“ aluminium to prevent the ball-bearing head from damage.

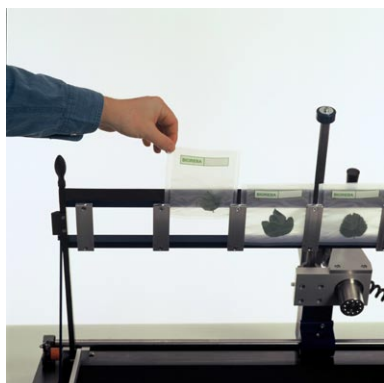
Store the HOMEX at a dry place to avoid corrosion. When not in use for several weeks/months, additionally wipe all metal parts with a soft cloth wetted with oil (or by using an anti corrosion protection spray).

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