HOMEX 7 USER GUIDE

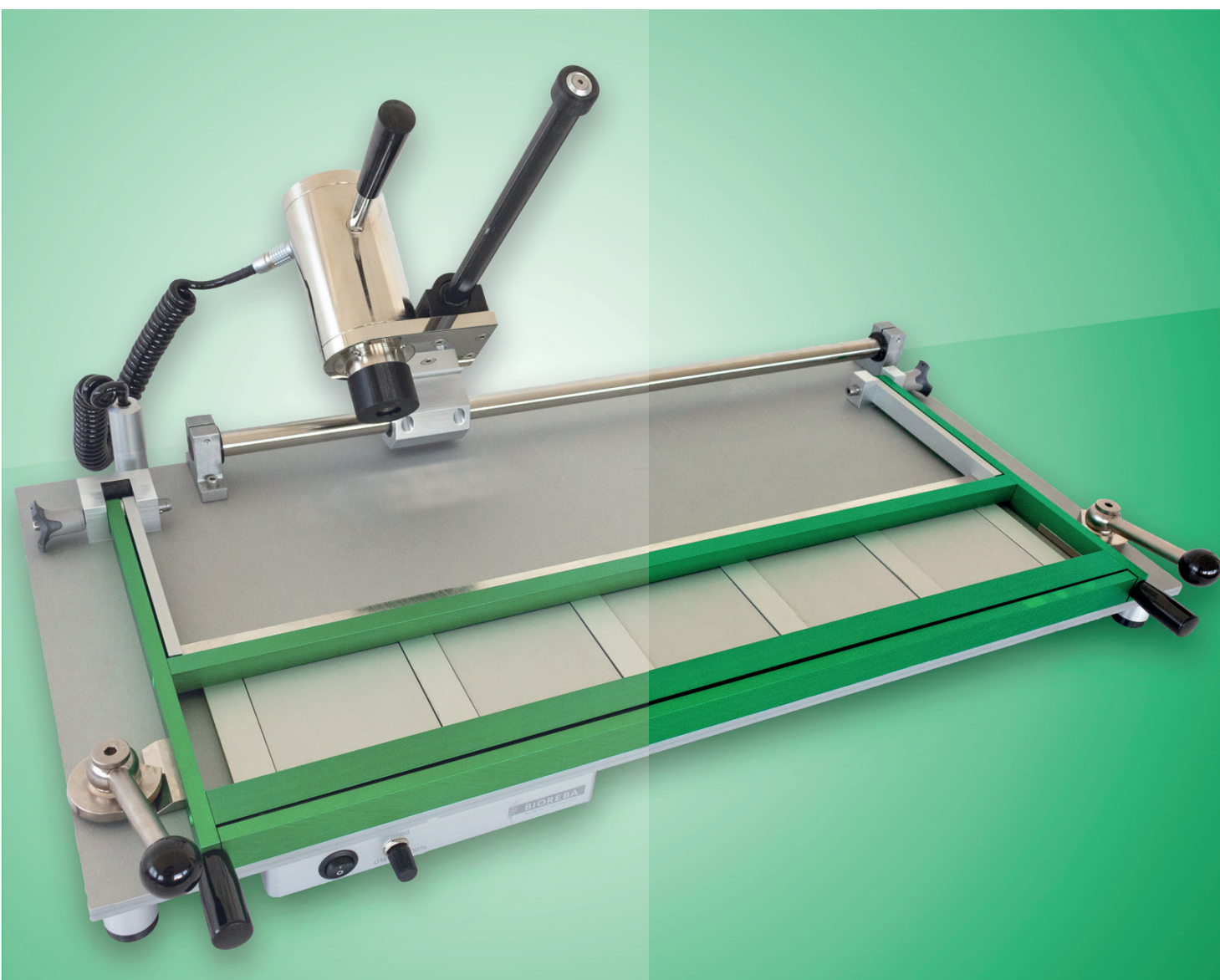
Table of contents

1	Foreword	3
1.1	Content check	3
2	Safety warning and guidelines	3
2.1	Safe Operating Procedures	3
3	Maintenance of the HOMEX 7	4
4	Warranty	5
5	Introduction	5
Table 1	Specifications	6
6	Operation guide	6
6.1	Replacing the rack	7
6.2	Replacing the ball-bearing head	7
6.3	RackClip: rack holder for more flexibility	8
7	Frequently Asked Questions	9
8	Packing list and Test certificate	9

Disposal

Observe the relevant legal regulations when disposing of the product.

Information on the disposal of electrical and electronic devices in the European Union:

Within the European Union, the disposal of electrical devices is regulated by national regulations based on EU Directive 2012/19/EU pertaining to waste electrical and electronic equipment (WEEE).

As the disposal regulations may differ from one country to another within the EU, please contact our local distributor for more information.

1. Foreword

Thank you for purchasing the HOMEX 7, a semi-automated homogenizing device for the grinding of plant samples such as leaves, wood, roots, seeds, and tubers.

The HOMEX 7 is ideal for high throughput applications and allows you to process up to 5 samples in rapid succession using ergonomic BIOREBA extraction bags. This user guide describes the HOMEX 7 functionalities, safety and operation instructions. Please read this user guide carefully before using the HOMEX 7 for the first time and keep it for later reference and troubleshooting.

1.1. Content check

Please check the packing list (see chapter 8) when you first unpack the HOMEX 7. If the instrument is damaged or items are missing, please contact us immediately.

2. Safety warnings and guidelines

Read this user guide before operating the instrument for the first time and follow the safe operating procedures described below.

2.1. Safe Operating Procedures

The operation, maintenance and repair of the instrument should comply with the basic usage guidelines and the warnings listed below. Failure to comply with the safe operating procedures may result in injury, affect the life span of the instrument and may result in the invalidation of the warranty.

- Do not carry the HOMEX 7 by yourself. Do not carry the HOMEX 7 holding it by the rack, the handle, or by any other external part. For short distances at least 2 people should carry the instrument holding it by the base plate.
 - If you need to ship the HOMEX 7, always use the original packaging.
 - Do not drop the HOMEX 7.
 - This device has been designed for indoor and non-extreme temperature conditions only.
 - In case the device is malfunctioning or is damaged, the operator should not try to repair it. This will result in the loss of warranty and may cause serious injury. If the device is not working properly or is defect, please contact the vendor immediately.
 - Before connecting the device to a power socket, make sure the voltage stated on the rating label of the product corresponds to the voltage of your power socket.
 - Ensure that the power plug and the power supply do not come into contact with water or moisture.
 - Do not disconnect the device from the power socket by pulling on the power cord. This could damage the power cord or the power socket. Disconnect the cord from the power socket by gently pulling the plug of the power cord.
 - Always place the instrument on a horizontal, stable and even surface. The surface must be resistant to heat and fluids.
 - Any liquid in contact with the electronic parts of the device may cause damage and provide a significant hazard.
 - The instrument should be placed in a dry, cool and dust-free environment with no exposure to direct sunlight. The instrument should be used only in a well-ventilated environment, with no corrosive gas or strong disturbing magnetic field, away from central heating and other heat sources.
 - When you finished to work switch off the instrument.
 - Pull out the power cord when you don't use the device for a long period of time. Clean the device and cover it with a cloth to protect from dust and sun.
- Do not:
- Use the instrument for anything other than its intended use.
 - Use a different power supply unit than the one supplied (Art. No. 400028).
 - Use the instrument if the cable to the power supply unit is damaged.
 - Contact the vendor in the following situations:
 - The instrument is drenched or fire burned.
 - The instrument produces abnormal noises or smells.
 - The instrument dropped.
 - The instrument does not function properly.

- If the power supply or the cable is damaged, it must be replaced by the manufacturer or companies approved by the manufacturer.

3. Maintenance of the HOMEX 7

The HOMEX 7 does not contain parts that rust. However, it is important to store and operate the HOMEX 7 in a dry place and to clean the instrument after each use. Prior to cleaning, power off the instrument and remove the connector plug. Immediately remove spills of extraction buffer with a humid soft cloth, then wipe dry - in particular when buffer has entered the ball-bearing head. Grease the two gliding axes (Fig. 1) after intense use or weekly with a soft cloth, wetted with oil or by using an anti-corrosion protection spray. Periodically check the holding-screws (especially those on the grinding head and on the rack) and tighten them if necessary. Replace the aluminium bars holding the bags when they are worn. They are made of "soft" aluminium to prevent the ball-bearing head from damage.

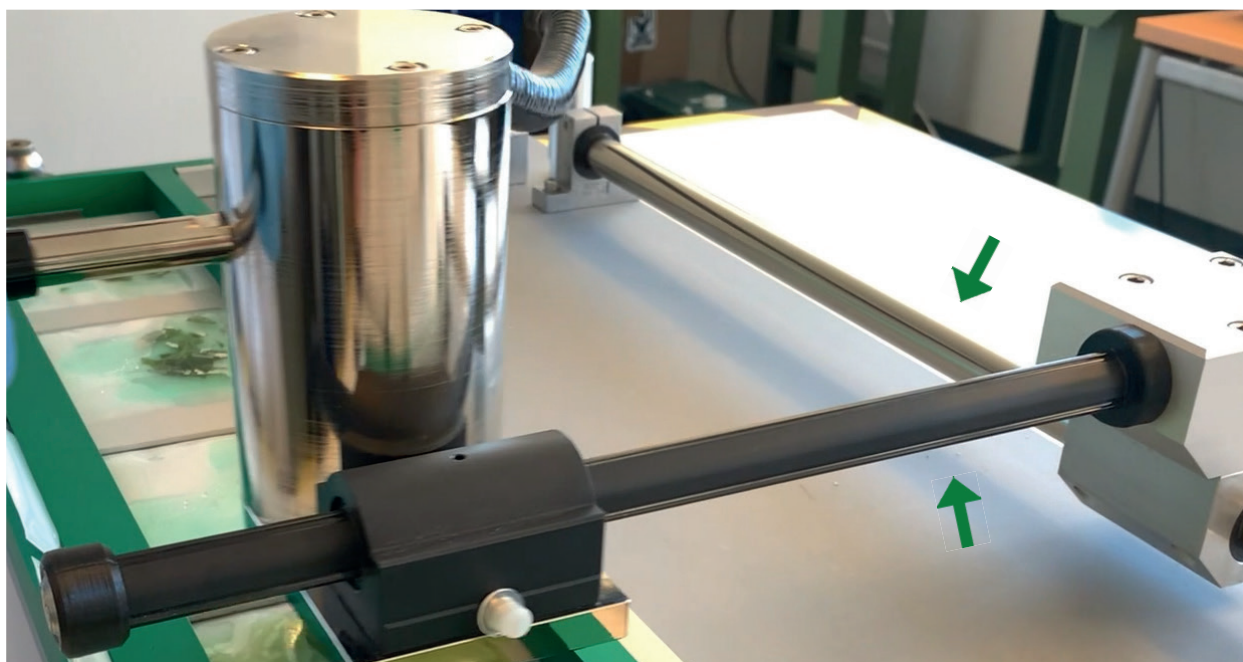
Fig. 1 Gliding axes

4. Warranty

Product Information

Warranty for BIOREBA equipment

BIOREBA equipment is of high quality and is checked for correct performance before shipment. The equipment is fully guaranteed against defective parts and materials, including defects caused by poor workmanship, for a period of 24 months from the date of delivery.

BIOREBA will repair or replace defective parts or materials during the term of warranty at no extra charge for materials and labor provided that the products were used and maintained in accordance with the manufacturer's instructions.

This guarantee is subject to the following exclusions:

- i. Any defects caused by normal wear and tear.
- ii. Defects caused by accident, misuse, misapplication, or as a result of incorrect power.
- iii. Defects caused by ecological disaster.

After delivery, please notify us (admin@bioreba.ch) or your distributor within 72 hours of receipt of the goods, of any damage, loss, or failure to meet product specifications, but also if a product fails to perform as expected during warranty, when used according to standard procedures. Please include with your notification the completed QS document "Complaint sheet for BIOREBA equipment" (form is available on our homepage www.bioreba.com).

All returns must be authorized prior to returning material.

Please return the equipment carefully packed - freight collect - to your distributor (or directly to Switzerland). BIOREBA AG will not be responsible for damage resulting from careless or insufficient packaging.

Products no longer under guarantee have to be sent back to Switzerland and will be repaired/serviced at your charge.

5. Introduction

The HOMEX 7 is a instrument designed for the efficient homogenization of plant tissues. Plant material including roots, leaves, and wood slices, is placed into a BIOREBA extraction bag with the appropriate amount of buffer. The standard rack can accomodate up to five BIOREBA "Universal" or "Universal-U-form" extraction bags, while the long rack can hold up to four „Universal-Long“ extraction bags for larger samples. The homogenized tissue is filtered through the synthetic intermediate layer of the BIOREBA bags, resulting in a particle-free homogenate that can be seamlessly used for downstream applications such as nucleic acid extraction, PCR, ELISA, Lateral Flow Immunochromatography, and more.

Although the HOMEX 7 homogenizes any kind of plant tissues, it may be necessary to pre-treat certain samples before homogenization: e.g., seeds may need prior soaking, while wood slices and thick/hard samples may need to be flattened with a rubber hammer (not provided).

The HOMEX 7 can process up to 250 samples per hour. The buffer volume should not exceed 4 ml for BIOREBA extraction bags "Universal-U-form" (Art. No. 440100), 10 ml for BIOREBA extraction bags "Universal" (Art. No. 430100), and 50 ml for BIOREBA extraction bags "Universal-Long" (Art. No. 470100).

Trademark	HOMEX
Model	HOMEX 7
Power supply unit	100 - 240 V AC, 50/60 Hz, 2A
Power	160 W
Rating	24 V, 6.67 A
Max speed	3000 rpm / adjustable (scale 1-10)
Sound level	70.5 dB (no sample) - 72.4 dB (with sample)
Dimensions	834 (L) x 451 (W) x 235 (H) mm Max height (rack or splined shaft in vertical position): 460 mm
Net weight (without power supply unit)	24.8 kg
Sample capacity	• 1-5 "Universal" or "Universal-U-form" using a Standard Rack (Art. No. 400014S) • 1-4 "Universal-Long" using a Long Rack (Art. No. 4000164)
BIOREBA Extraction bags	• "Universal" (Art. No. 430100) • max sample volume: 1 g • max buffer volume: 10 ml • dimensions: 120 x 150 mm • "Universal-Long" (Art. No. 470100) • max sample volume: 5 g • max buffer volume: 50 ml • dimensions: 150 x 280 mm • "Universal-U-form" (Art. No. 440100) • designed for seeds • max sample volume: 0.5 g • max buffer volume: 4 ml • dimensions: 120 x 150 mm

Table 1. Specifications

6. Operation guide

Follow these steps to homogenize up to 5 samples or watch a video tutorial by scanning following QR code:

Scan me

- First connect the power unit to the HOMEX 7 and then plug the power unit into the electrical outlet. Make sure that all cables are plugged in tightly.
- Place the plant material into BIORÉBA extraction bags along with the appropriate amount of buffer. If necessary, pre-treat samples such as seeds and wood slices before homogenizing.
- Insert up to 5 extraction bags into the rack (Fig. 2).
- Close the rack and pull the two tension levers (Fig. 3) towards the front of the instrument to securely lock the extraction bags in the homogenizing position.
- Set the speed from 0 to 100% using the potentiometer (Fig. 4).
- Press the "I/O" button (Fig. 4) to begin homogenizing the samples. Glide the grinding head over the bags (Fig. 5) without applying excessive pressure. Ensure you keep a firm grip on the handle to prevent the homogenizing apparatus from slipping.

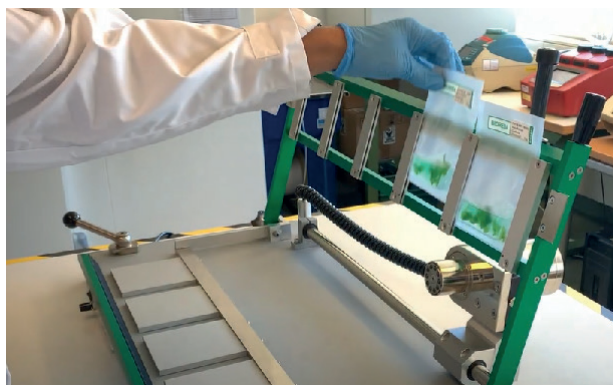
Fig. 2

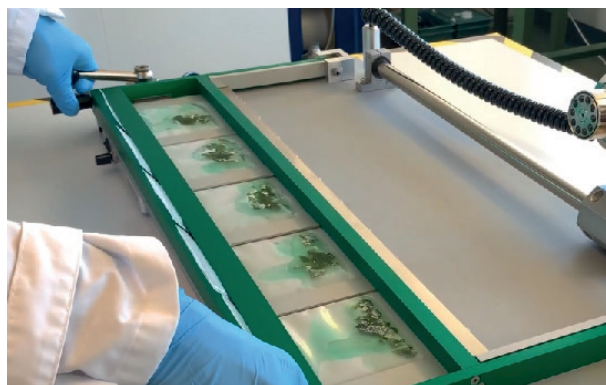
Fig. 3

Fig. 4

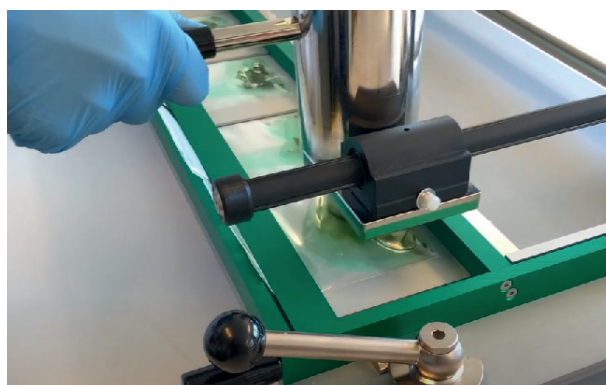
Fig. 5

6.1 Replacing the rack

To replace the rack, press the knobs at the two fixing screws and pull the screws out (Fig. 6). Insert the new rack and reposition the screws. When moving or carrying the rack, exercise caution, as the folding parts may clamp your fingers. Watch a video tutorial on how to change the rack by scanning following QR code.

Video: How to
change the rack
for HOMEX 7

Fig. 6

Video: How to
remove the ball-
bearing head for
HOMEX 7

6.2 Replacing the ball-bearing head

To replace or clean the ball-bearing head, first use the provided 2 mm Allen key to remove the four screws on the side, and then remove the cover case (Fig. 7). To replace the ball-bearing head entirely, use the 2 mm Allen key to pull out the head fixation screw (Fig. 8). Insert the new ball-bearing head. Apply Loctite 243 (threadlocker, not included) to the fixation screw (Fig. 9) and insert it towards the flat surface of the metallic shaft, as shown in Fig 10. Loctite will be effective after 10 minutes. However, it is preferable to wait 24 hours before using the homogenizer again. Watch a video tutorial on how to remove and replace the ball-bearing head by scanning the QR code above.

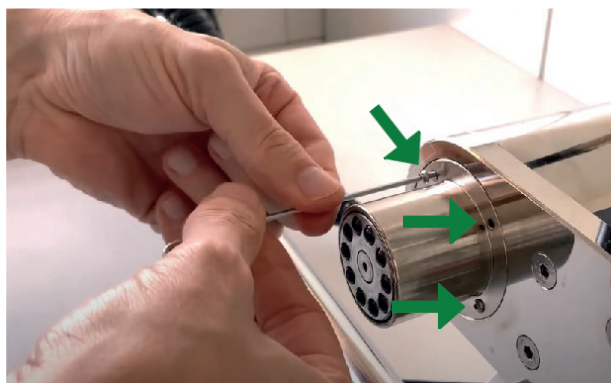
Fig. 7

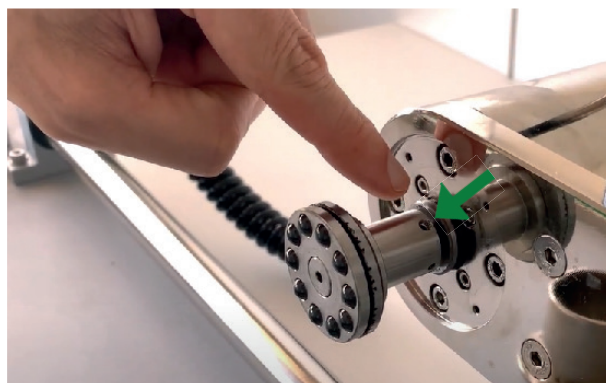
Fig. 8

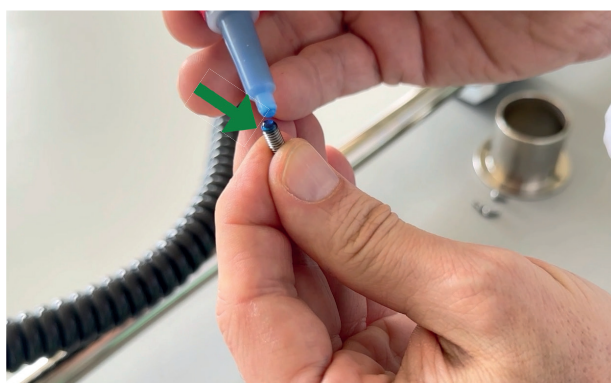
Fig. 9

Fig. 10

6.3 RackClip: rack holder for more flexibility

The RackClip (rack holder) simplifies the insertion of BIOREBA extraction bags into the open standard rack. Mount the RackClip onto the rack using a 2.5 mm Allen key. The tilt angle of the rack is determined by the position of the RackClip. When the RackClip is open (Fig. 11, top), you can open and close the rack freely. When the RackClip is closed (Fig. 11, bottom), the rack remains tilted, facilitating the insertion of extraction bags (Fig. 12). Watch the video tutorial on how to assemble and use the RackClip by scanning the following QR code:

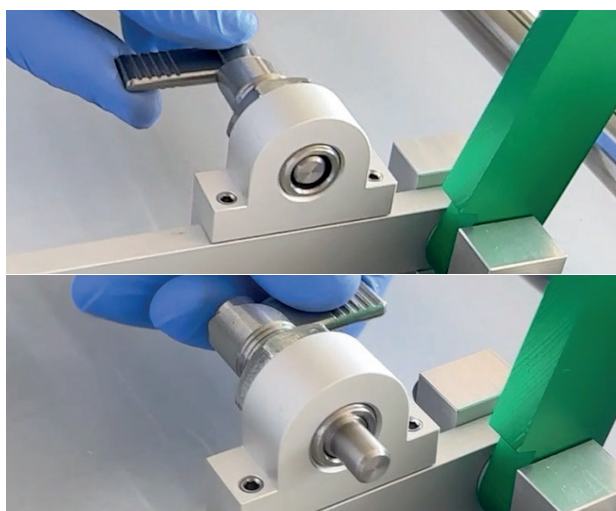
Fig. 11

Fig. 12

Scan me

Video: HOMEX RackClip:
installation and use

7. Frequently Asked Questions

Problem / Question	Possible cause	Solution
The instrument won't start after pressing the I/O button.	- The cables are not connected properly. - The cables or the power supply unit are damaged. - Mechanic parts of the instrument may be defective or damaged.	- When you plug the cable into the HOMEX 7 socket, you should hear a „click“ - Replace the cables or power supply unit with one provided by BIOREBA. - Do not further use the instrument and contact admin@bioreba.ch.
Liquid is spilling out of the extraction bag during homogenization.	- Buffer and/or sample volume has been exceeded. - The rack has not been closed tightly (Fig. 3).	- Do not exceed the volumes described in table 1. - Pull the tension levers tightly toward the front of the HOMEX 7 (Fig. 3).
The extraction bag breaks during homogenization.	- Too much pressure has been applied to the sample. - The sample may be too hard/edgy (e.g., wood samples, hard roots or big seeds) causing the bag to break. - Mechanic parts of the instrument (e.g. ball-bearing head) or the extraction bags may be defect or damaged.	- Avoid applying excessive pressure to the sample. The weight of the grinding head should be sufficient to homogenize most samples. - Pre-treat hard samples with a rubber hammer or pre-soak them in buffer for several hours before homogenization. - Consider replacing the ball-bearing head if you notice any defect.
The sample is not homogenized appropriately.	Some samples such as seeds or woody samples may need pre-treatment.	- For seeds it is recommended to soak them in buffer for at least 4 hours prior to homogenization. - For woody samples a pre-treatment with a rubber hammer is recommended. - Try to modify the speed of homogenization to obtain better results.

8. Packing list and replacement parts

Standard rack version

- 1 x HOMEX 7 incl. Standard Rack (Art. No 400007S)
- 1 x RackClip (rack holder) (Art. No. 400037)
- 100 x BIOREBA extraction bags "Universal" (Art. No. 430100)
- 100 x BIOREBA extraction bags "Universal-U-form" (Art. No. 440100)

Long rack version

- 1 x HOMEX 7 incl. Long Rack (Art. No. 4000074)
- 100 x BIOREBA extraction bags "Universal-Long" (Art. No. 470100)

Both versions also contain

- 1 x Power supply unit (Art. No. 400028)
- 1 x Cable to the power supply unit (Art. No. 400038)
- 1 x Allen key 2 mm (Art. No 545831)
- 1 x Allen key 2,5 mm (Art. No 545832)
- 1 x Oil bottle (Art. No 450160)
- 1 x User Guide

Replacement parts (not included in the packing list)

- Standard Rack for HOMEX 7 (Art. No. 400014S)
- Long Rack for HOMEX 7 (Art. No. 4000164)
- Replacement head for HOMEX 7 (Art. No. 400017)
- Aluminium bars (6x) for Standard Rack HOMEX 7 (Art. No. 400027)

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**To which this declaration relates is in conformity with the following standard(s) or other
normative document(s).**

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IEC 61010-1:2010 + AMD1:2016
EN 61010-1:2010 + A1:2019
IEC 61010-2-051:2015

EN 61010-2-051:2015
EU Group Differences, US & CA National Differences

(Titel und/oder Nummer sowie Ausgabedatum der Norm(en) oder der anderen normativen Dokumente)
(Title and/or number and date of issue of the standard(s) or other normative document(s))
Titre et ou no et date de publication de la (des) norme(s) ou autre(s) document(s) normatif(s)

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Following the provisions of Directive(s) (if applicable):

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