HOMEX 7 USER GUIDE

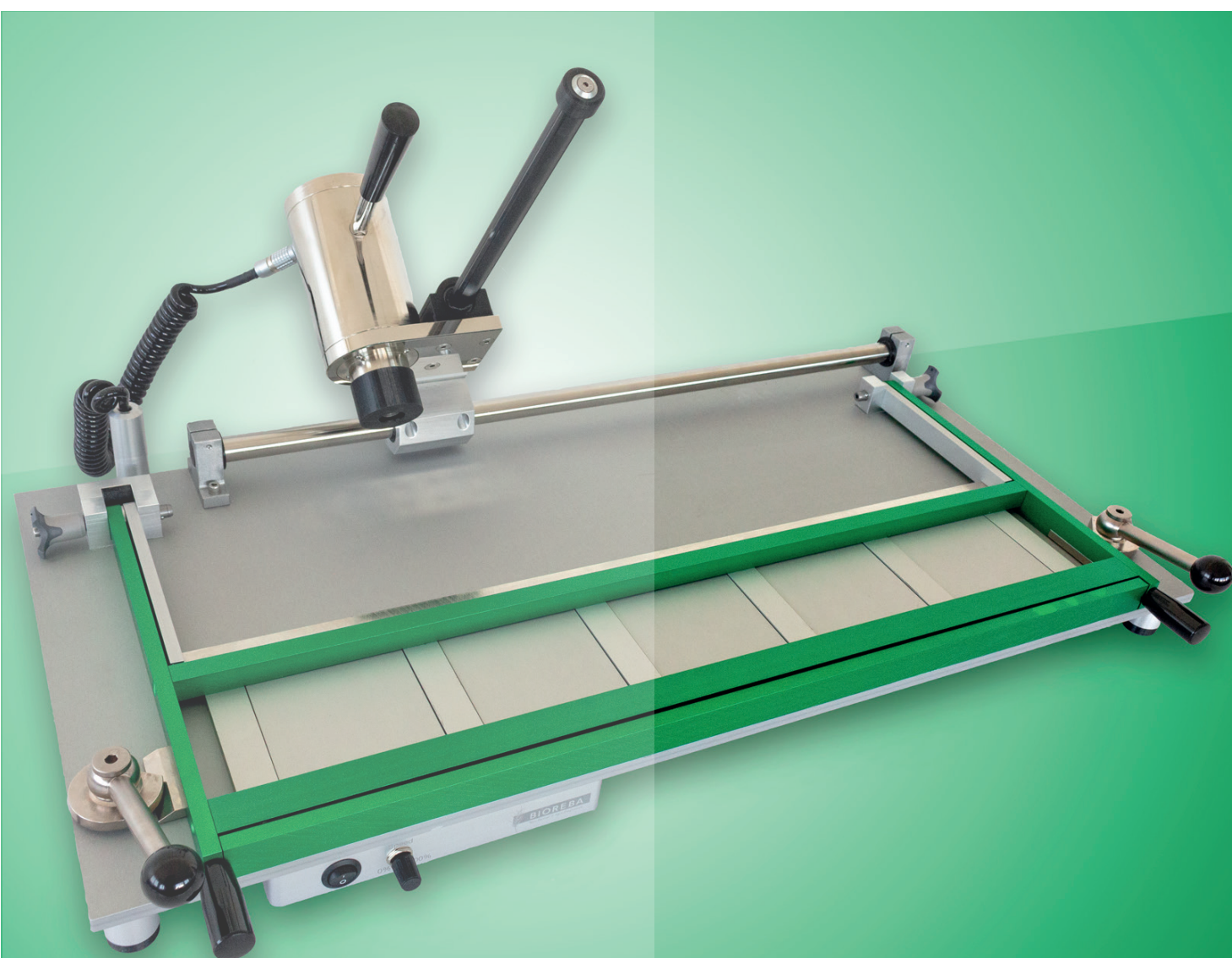
Table of contents

1	Foreword	3
1.1	Content check	3
2	Safety warning and guidelines	3
2.1	Safe Operating Procedures	3
3	Maintenance of the instrument	4
4	Warranty	5
5	Introduction	5
Table 1.	Specifications	6
6	Operation guide	6
5.1	Replacing the rack	7
5.2	Removing the ball-bearing head	7
7	Frequently Asked Questions	8
8	Packing list and Test certificate	9

1. Foreword

Thank you for purchasing the HOMEX 7, a semi-automated homogenizing device for the grinding of plant samples such as leaves, wood, roots, seeds, and tubers.

The HOMEX 7 is ideal for high throughput applications and allows you to load up to 5 extraction bags. This user guide describes the HOMEX 7 functionalities, safety and operation instructions. Please read this user guide carefully before using the HOMEX 7 for the first time and keep it for later reference and troubleshooting.

1.1. Content check

Please check the packing list (see chapter 7) when you first open the HOMEX 7 packaging. If the instrument is damaged or items are missing, please contact us immediately.

2. Safety warnings and guidelines

Read this user guide before operating the instrument for the first time and follow the safe operating procedures described below.

2.1. Safe Operating Procedures

The operation, maintenance and repair of the instrument should comply with the basic usage guidelines and the warnings listed below. Failure to comply with these guidelines may result in injury, affect the life span of the instrument and may result in the invalidation of the warranty.

- Do not carry the HOMEX 7 by yourself. Never carry the HOMEX 7 holding it by the rack, the handle, or by any other external part. For short distances at least 2 people should carry the instrument holding it by the base plate.
 - In case you need to ship the HOMEX 7, always use the original packaging.
 - Do not drop the HOMEX 7.
 - This device has been designed for indoor and non-extreme temperature conditions use only.
 - In case the device is malfunctioning or is damaged, the operator should not try to repair it. This will result in the loss of warranty and may cause serious injury. If the device is not working properly or is defective, please contact the manufacturer immediately.
 - Before connecting the device to a power socket, make sure the voltage stated on the rating label of the product corresponds to the voltage in your power socket.
 - Ensure that the power plug does not come into contact with water or moisture.
 - Do not disconnect the device from the power socket by pulling on the power cord. This could damage the power cord or the power socket. Disconnect the cord from the power socket by gently pulling the plug of the power cord.
 - Always place the instrument on a horizontal, stable and even surface. The surface must be resistant to heat and fluids.
 - Any liquid in contact with the electronic parts of the device may cause damage and provide a significant hazard.
 - The instrument should be placed in a dry, cool and dust-free environment with no exposure to direct sunlight. The instrument should be used only in a well-ventilated environment, with no corrosive gas or strong disturbing magnetic field, away from central heating and other heat sources.
 - Switch off the instrument when you finished to work.
 - Pull out the connector plug when you don't use the device for a long period of time. Clean the instrument and cover it with a cloth to protect from dust.
- Do not:
- Use the instrument for anything other than its intended use.
 - Use a different power supply unit than the one supplied (Art. No. 400028).
 - Use the instrument if the cable to the power supply unit is damaged.
 - Pull the connector plug and contact the vendor in the case:
 - The instrument is drenched or fire burned.
 - The instrument produces abnormal noises or smells.
 - The instrument dropped.
 - The instrument does not function properly.

- Keep the connector plug away from heat and damp. If the supply cord is damaged, it must be replaced by the manufacturer or companies approved by the manufacturer.

3. Maintenance of the instrument

The HOMEX 7 does not contain parts which may rust. However, it is important to store and operate the HOMEX 7 at a dry place and clean the instrument after use. Prior to cleaning, power off the instrument and remove the connector plug. Immediately remove spills of extraction buffer with a humid soft cloth, then wipe dry - in particular when buffer entered the ball-bearing head. Grease the two gliding axes (Fig. 1) daily with a soft cloth, wetted with oil or by using an anti-corrosion protection spray. Check the holding-screws of the motor and the rack and tighten them if necessary. Replace the aluminium bars holding the bags when they look worn. They are made of "soft" aluminium to prevent the ball-bearing head from damage.

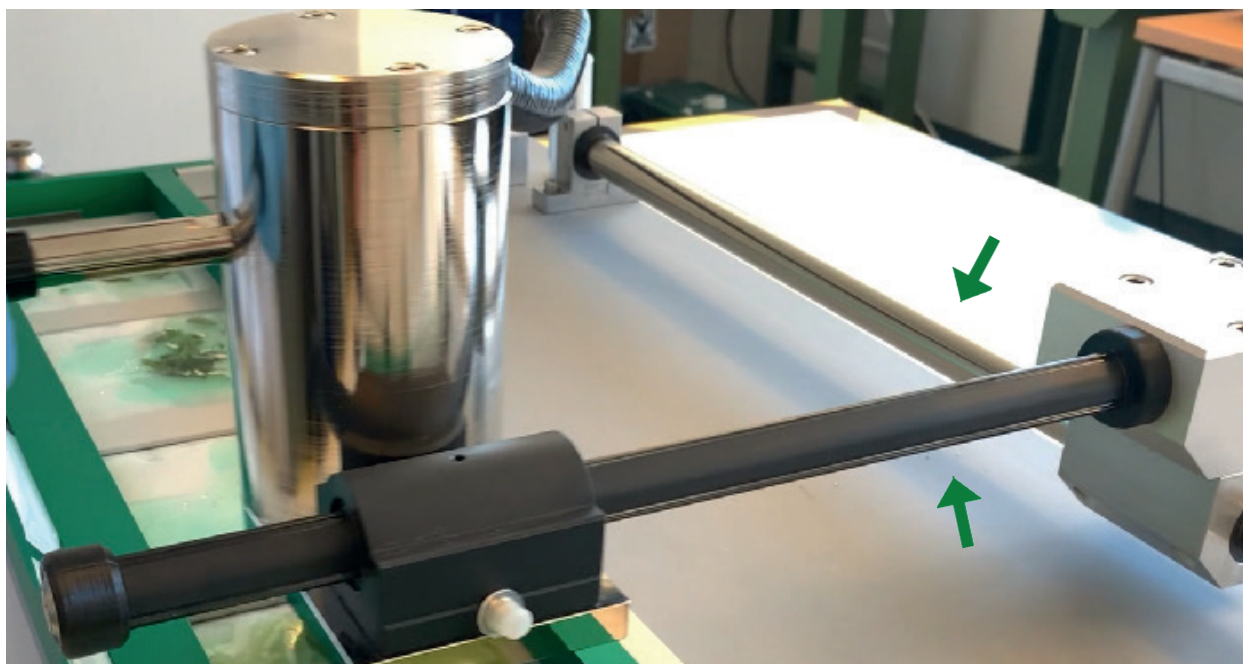
Fig. 1 Gliding axes

4. Warranty

Product Information

Warranty for BIOREBA equipment

BIOREBA equipment is of high quality and is checked for correct performance before shipment. The equipment is fully guaranteed against defective parts and materials, including defects caused by poor workmanship, for a period of 24 months from date of delivery.

BIOREBA will repair or replace defective parts or materials during the term of warranty at no extra charge for materials and labor provided that the products were used and maintained in accordance with the manufacturer's instructions.

This guarantee is subject to the following exclusions:

- i. Any defects caused by normal wear and tear.
- ii. Defects caused by accident, misuse, misapplication or as result of incorrect power.
- iii. Defects caused by ecological disaster.

After delivery, please notify us (admin@bioreba.ch) or your distributor within 72 hours of receipt of the goods, of any damage, loss or failure to meet product specifications, but also if a product fails to perform as expected during warranty, when used according to standard procedures. Please include with your notification the completed QS document "Complaint sheet for BIOREBA equipment" (form is available on our homepage www.bioreba.com).

All returns must be authorized prior to returning material.

Please return the equipment carefully packed - freight collect - to your distributor (or directly to Switzerland). BIOREBA AG will not be responsible for damage resulting from careless or insufficient packaging.

Products no longer under guarantee have to be sent back to Switzerland and will be repaired/serviced at your charge.

5. Introduction

The HOMEX 7 is an instrument for the homogenization of plant tissues. Plant material such as roots, leaves, and wood slices, is inserted into a BIOREBA extraction bag with the appropriate amount of buffer. Up to 5 BIOREBA extraction bags "Universal" or "Universal-U-form" can be inserted into the standard rack. For bigger samples, up to 4 BIOREBA extraction bags "Universal-Long" can be inserted into a 4/25 rack. The homogenized tissue is filtered by the synthetic intermediate layer of the BIOREBA bags and can then be easily used for downstream applications such as nucleic acid extraction, PCR, ELISA, Lateral Flow Immunochromatography, etc.

Although the HOMEX 7 homogenizes any kind of plant tissues, it may be necessary to pre-treat certain samples before homogenization: e.g., seeds may need presoaking, while wood slices and thick samples may need to be flattened with a rubber hammer (not provided).

The HOMEX 7 can process up to 250 samples per hour. The buffer volume should not exceed 4 ml for BIOREBA extraction bags "Universal-U-form" (Art. No. 440100), 10 ml for BIOREBA extraction bags "Universal" (Art. No. 430100), and 50 ml for BIOREBA extraction bags "Universal-Long" (Art. No. 470100).

Trademark	HOMEX
Model	HOMEX 7
Power supply unit	100 - 240 V AC, 50/60 Hz, 2A
Power	160 W
Rating	24 V, 6.67 A
Max speed	3000 rpm / adjustable (scale 1-10)
Sound level	70.5 dB (no sample) - 72.4 dB (with sample)
Dimensions	834 x 451 x 235 mm
Weight (without power supply unit)	24.8 kg
Sample capacity	• 1-5 using a Standard rack (Art. No. 400014S) - or • 1-4 using a Rack 4/25 (Art. No. 4000164)
BIOREBA Extraction bags	• "Universal" (Art. No. 430100) • max sample volume: 1 g • max buffer volume: 10 ml • dimensions: 120 x 150 mm • "Universal-Long" (Art. No. 470100) • max sample volume: 5 g • max buffer volume: 50 ml • dimensions: 150 x 280 mm • "Universal-U-form" (Art. No. 440100) • designed for seeds • max sample volume: 0.5 g • max buffer volume: 4 ml • dimensions: 120 x 150 mm

Table 1. Specifications

6. Operation guide

Follow these steps to homogenize up to 5 samples or watch a video tutorial by scanning following QR code:

- Plug the power supply unit into a socket and connect the HOMEX 7 with the power supply cable at the front left side of the instrument.
- Prepare the samples using the BIOREBA extraction bags and add the appropriate amount of buffer. If necessary, pre-treat samples such as seeds and wood slices before homogenizing.
- Insert up to 5 extraction bags in the open rack (see Fig. 2).
- Close the rack and pull the two tension levers (see Fig. 3) toward the instrument's front side in order to lock the extraction bags in the homogenizing position.
- Set the speed using the potentiometer (see Fig. 4).
- Press the "I/O" button (Fig. 4) and start homogenizing the samples by gliding the "ball-bearing head" over the bags (see Fig. 5) without applying too much pressure. Keep the handle firmly in your hand in order not to drop the homogenizing apparatus.

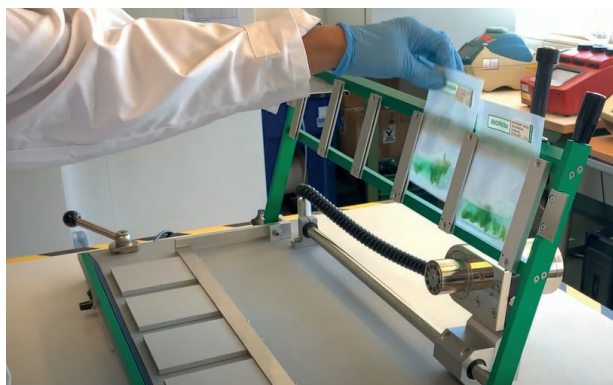
Fig. 2

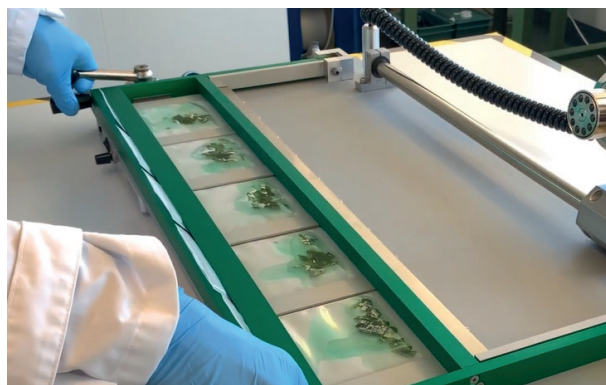
Fig. 3

Fig. 4

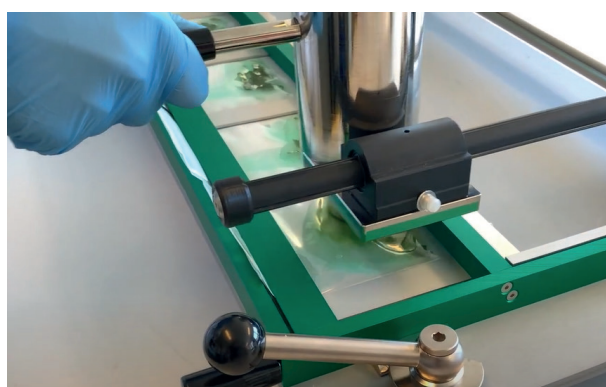
Fig. 5

6.1 Replacing the rack

In order to replace the rack, you need to push the knobs (Fig. 6) at the two fixing screws and pull the screws out. Insert the new rack and put the screws back in position. When moving or carrying the rack, try to be careful as the folding parts may clamp your fingers. Watch a video tutorial on how to change the rack by scanning following QR code.

Video: How to change the rack for HOMEX 7

Fig. 6

Video: How to remove the ball-bearing head for HOMEX 7

6.2 Removing the ball-bearing head

In order to replace or clean the ball-bearing head, you first need to remove the 4 screws on the side using the provided 2 mm allen key and then remove the cover-case (Fig. 7). To remove the ball-bearing head entirely, the head fixation screw has to be pulled out as well using a 2 mm allen key (Fig. 8). Watch a video tutorial on how to change the ball-bearing head by scanning QR code above.

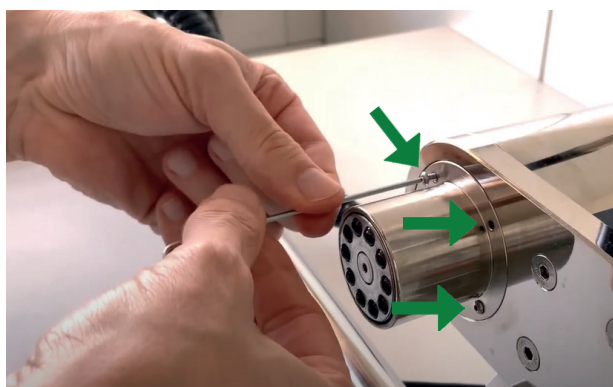
Fig. 7

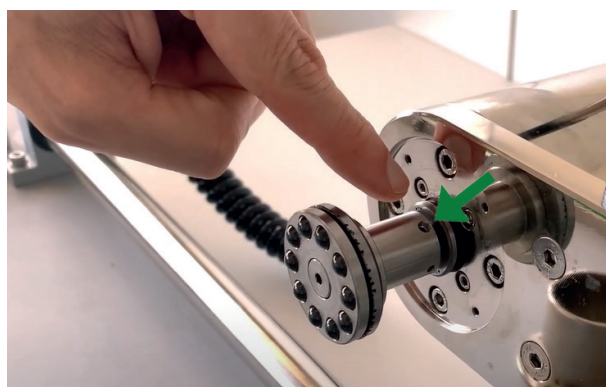
Fig. 8

7. Frequently Asked Questions

Problem / Question	Possible cause	Solution
Instrument won't start after pressing the I/O button.	• Cable is not connected properly. • Cable or power supply unit are damaged. • Mechanic parts of the instrument may be defect or damaged.	• Replace cable or power supply unit with one provided by BIOREBA. • Do not further use the instrument and contact admin@bioreba.ch.
Homogenate is spilling out of the extraction bag during homogenization.	• Buffer and/or sample volume has been exceeded. • The rack has not been closed tightly (Fig. 3).	• Do not exceed the volumes described in table 1. • Pull the clamping-screws tightly (Fig. 3).
Extraction bag breaks during homogenization.	• Too much pressure has been applied to the sample. • Sample may be too hard/edgy (e.g., wood samples, hard roots or big seeds).	• Avoid applying pressure to the sample. The weight of the homogenizing head should be enough to homogenize most of the samples. • Pre-treat hard samples with a rubber hammer or pre-soak them in buffer for several hours before homogenization.
Sample is not homogenized appropriately.	• Some samples such as seeds or woody samples may need pre-treatment.	• For seeds it is recommended to soak them in buffer for at least 4 hours prior to homogenization. • For woody samples a pre-treatment with a rubber hammer is recommended. • Try to modify the speed of homogenization to obtain better results.

8. Packing list

Standard rack version

- 1x HOMEX 7 incl. Standard Rack (Art. No 400007S)
- 100x BIOREBA extraction bags "Universal" (Art. No. 430100)
- 100x BIOREBA extraction bags "Universal-U-form" (Art. No. 440100)

4/25 rack version


- 1x HOMEX 7 incl. Rack 4/25 (Art. No. 4000074)
- 100x BIOREBA extraction bags "Universal-Long" (Art. No. 470100)

Both versions also contain

- 1x Power supply unit (Art. No. 400028)
- 1x Cable to the power supply unit (Art. No. 400038)
- 1x Allen key 2 mm (Art. No 545831)
- 1x Allen key 2,5 mm (Art. No 545832) (for the aluminum bars)
- 1x Oil bottle (Art. No 450160)
- 1x User Guide

 emc testcenter.zürich.ag				
EMC-Testcenter AG Moosackerstrasse 77 8105 Regensdorf SWITZERLAND Phone +41 44 302 45 00 E-mail info@emc-testcenter.com Homepage www.emc-testcenter.com	Accredited according to ISO / IEC 17025 Swiss Accreditation Service SAS (STS 0034) IECEE CBTL (TL0651) Listed according to Article 17 TGV at ASTRA  			
TEST CERTIFICATE EMCKC4170B				
Product :	Tissue homogenizer			
Trade mark :	Homex			
Type :	HOMEX 7			
Test results :	EMCKP4170.1B EMCKP4170.1B.ND EMCKP4170.2B			
A sample of this product was tested and found to be in conformity according to the following standards: IEC 61010-1:2010 + AMD1:2016 EN 61010-1:2010 + A1:2019 IEC 61010-2-051:2015 EN 61010-2-051:2015 EU Group Differences, US & CA National Differences				
Regensdorf, 2022-05-24	EMC-TESTCENTER AG  Roger MARTI Head of SAFETY Lab			
<table border="1" style="width: 100%; border-collapse: collapse;"> <tr> <td style="width: 33%; padding: 2px;">Date of Issue: 2022-05-24</td> <td style="width: 33%; padding: 2px;">Certificate No.: EMCKC4170B</td> <td style="width: 33%; padding: 2px;">Page 1 of 1</td> </tr> </table>		Date of Issue: 2022-05-24	Certificate No.: EMCKC4170B	Page 1 of 1
Date of Issue: 2022-05-24	Certificate No.: EMCKC4170B	Page 1 of 1		

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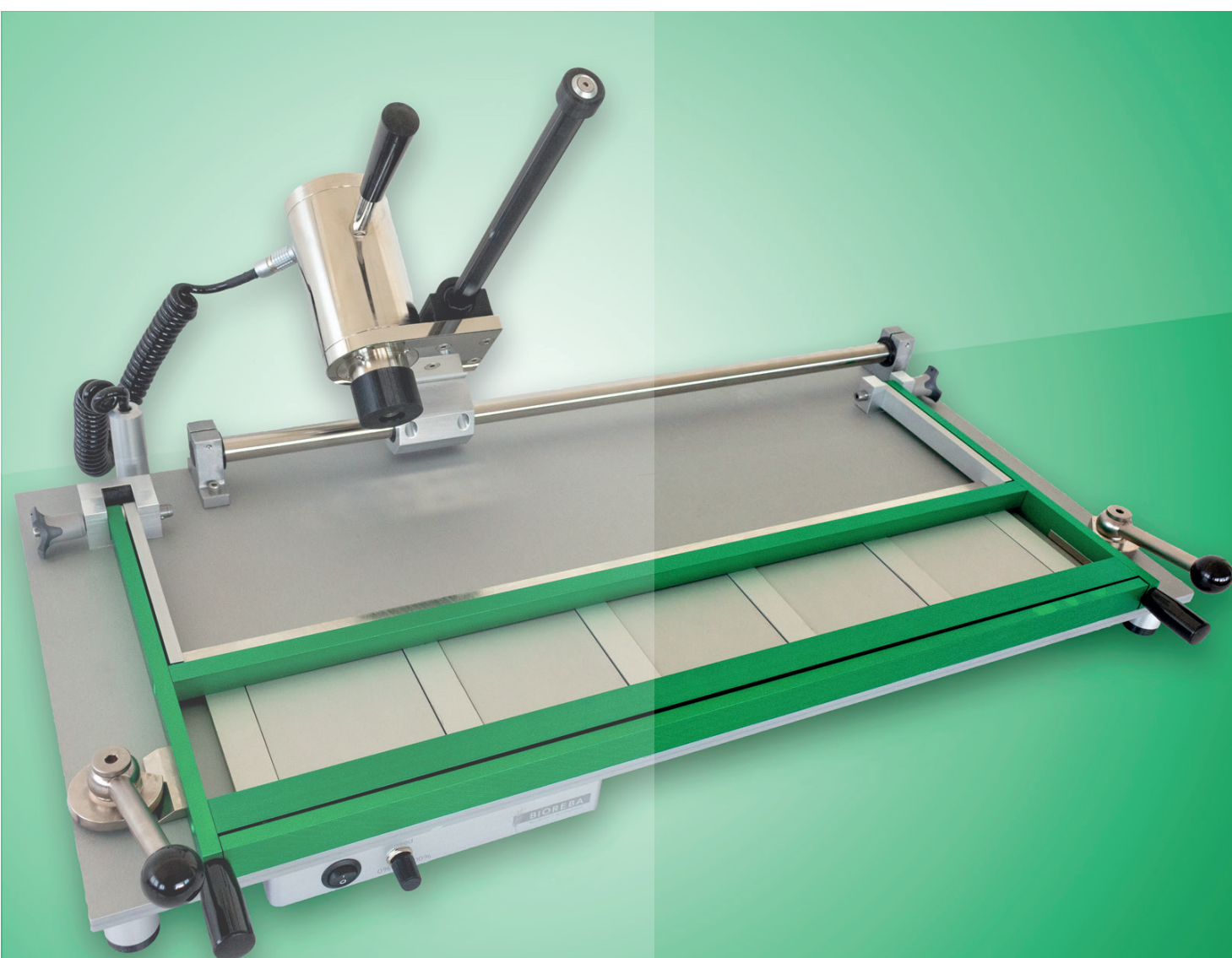 emc testcenter.zürich.ag				
EMC-Testcenter AG Moosackerstrasse 77 8105 Regensdorf SWITZERLAND Phone +41 44 302 45 00 E-mail info@emc-testcenter.com Homepage www.emc-testcenter.com	Accredited according to ISO / IEC 17025 Swiss Accreditation Service SAS (STS 0034) IECEE CBTL (TL0651) Listed according to Article 17 TGV at ASTRA  			
TEST CERTIFICATE EMCKC4170B				
Product :	Tissue homogenizer			
Trade mark :	Homex			
Type :	HOMEX 7			
Test results :	EMCKP4170.1B EMCKP4170.1B.ND EMCKP4170.2B			
A sample of this product was tested and found to be in conformity according to the following standards: IEC 61010-1:2010 + AMD1:2016 EN 61010-1:2010 + A1:2019 IEC 61010-2-051:2015 EN 61010-2-051:2015 EU Group Differences, US & CA National Differences				
Regensdorf, 2022-05-24	EMC-TESTCENTER AG  Roger MARTI Head of SAFETY Lab			
<table border="1" style="width: 100%; border-collapse: collapse;"> <tr> <td style="width: 33%;">Date of Issue: 2022-05-24</td> <td style="width: 33%;">Certificate No.: EMCKC4170B</td> <td style="width: 33%;">Page 1 of 1</td> </tr> </table>		Date of Issue: 2022-05-24	Certificate No.: EMCKC4170B	Page 1 of 1
Date of Issue: 2022-05-24	Certificate No.: EMCKC4170B	Page 1 of 1		

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emc testcenter.zürich.ag

EMC-Testcenter AG
Moosackerstrasse 77
8105 Regensdorf
SWITZERLAND

Phone +41 44 302 45 00
E-mail info@emc-testcenter.com
Homepage www.emc-testcenter.com

Accredited according to ISO / IEC 17025
Swiss Accreditation Service SAS (STS 0034)
IECEE CBTL (TL0651)
Listed according to Article 17 TGV at ASTRA

TEST CERTIFICATE EMCKC4170B

Product : Tissue homogenizer

Trade mark : Homex

Type : HOMEX 7

Test results : EMCKP4170.1B
EMCKP4170.1B.ND
EMCKP4170.2B

A sample of this product was tested and found to be in conformity according to the following standards:

IEC 61010-1:2010 + AMD1:2016
EN 61010-1:2010 + A1:2019
IEC 61010-2-051:2015
EN 61010-2-051:2015
EU Group Differences, US & CA National Differences

Regensdorf, 2022-05-24

EMC-TESTCENTER AG

Roger MARTI
Head of SAFETY Lab

Date of Issue: 2022-05-24	Certificate No.: EMCKC4170B	Page 1 of 1
---------------------------	-----------------------------	-------------

Your Partner in Agro-Diagnostics

www.bioreba.com