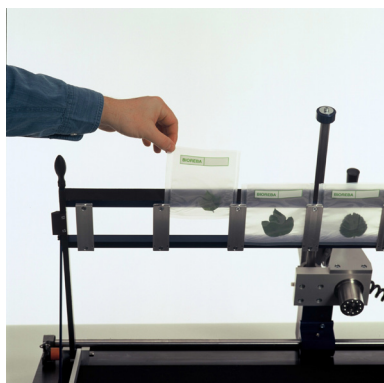
Product Information

HOMEX 6 Instructions

Filling the rack

Insert 1 to 5 extraction bags in the open rack.

Closing the rack

Close the rack by turning the two clamping-screws (1) to fix the bags in the homogenizing position.

Homogenizing

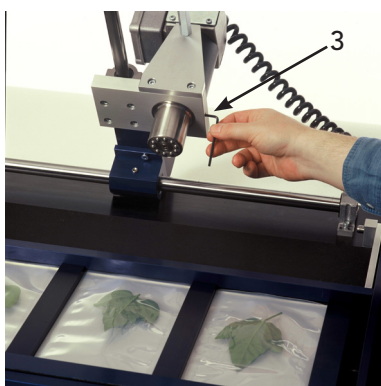
Switch the motor on and homogenize the sample by gliding the „ball-bearing head“ over the bags, using the handle.

Changing the rack

To change the rack, push the knobs (2) at the two fixing-screws and pull the screws out. Insert the rack and put the screws back in position.


Removing the ball-bearing head

To replace or clean the ball head, first unscrew the side screw (3) and then remove the cover. To completely remove the head, remove the head locking screw (4). When reassembling the head, it is important to insert the fixing screw (4) in the recess of the metal support. Scan the QR code to watch a short video tutorial (starts at 0:49).

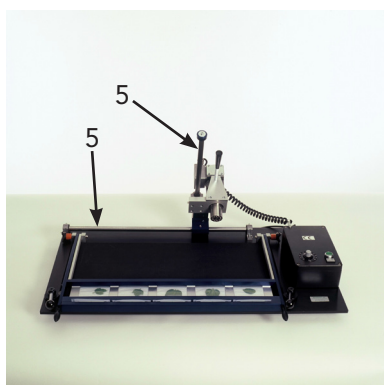Maintenance, storage and corrosion protection

Scan me.

Watch the tutorial on our YouTube channel.

Immediately remove spills of extraction buffer with a humid soft cloth, then wipe dry - in particular when buffer entered the ball-bearing head. Lubrify the two gliding axes (5) daily with a soft cloth, wetted with oil or by using an anti-corrosion protection spray. Check the holding-screws of the motor and the rack and tighten them if necessary. Replace the aluminium bars holding the bags when they look worn. They are made of „soft“ aluminium to prevent the ball-bearing head from damage.

HOMEX 6 contains parts which may rust. Store and operate the HOMEX 6 at a dry place to avoid corrosion.

1/1

Version: 5 - 05.02.2024

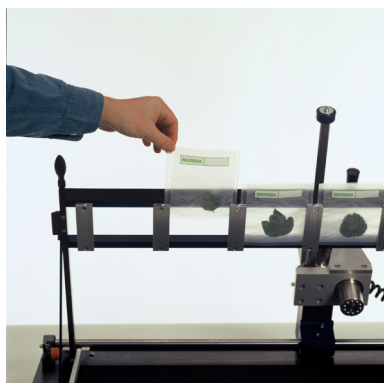
Adaptations from last version: Minor change in maintenance, new QR code for video tutorial.

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