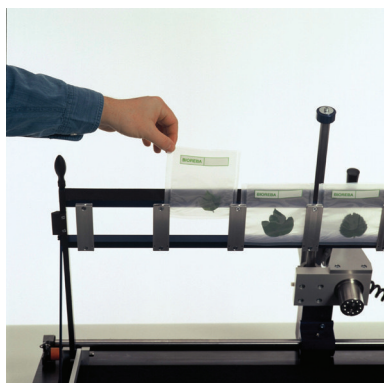
Product Information

HOMEX 6 Instructions

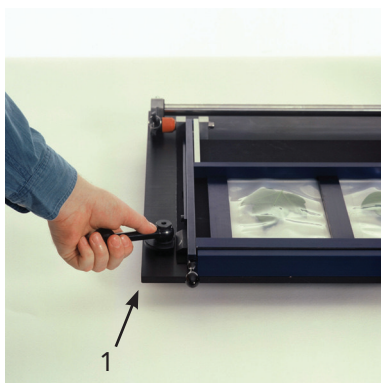
Filling the rack

Insert 1 to 5 extraction bags in the open rack.

Closing the rack

Close the rack by turning the two clamping-screws (1) to fix the bags in the homogenizing position.

Homogenizing

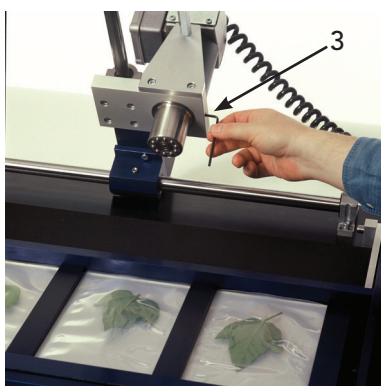
Switch the motor on and homogenize the sample by gliding the „ball-bearing head“ over the bags, using the handle.

Changing the rack

To change the rack, push the knobs (2) at the two fixing-screws and pull the screws out. Insert the rack and put the screws back in position.


Removing the ball-bearing head

To change or clean the ball-bearing head, first open the screw on the side (3) then remove the cover-case. To remove the ball-bearing head entirely, the head fixation screw (4) has to be pulled out as well.

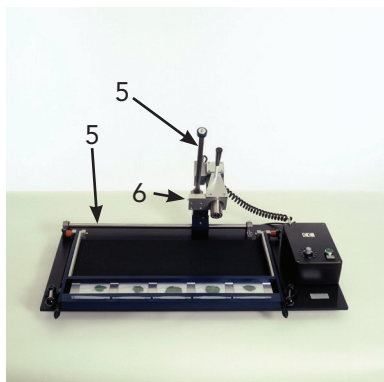Maintenance, storage and corrosion protection

Immediately remove spills of extraction buffer with a humid soft cloth, then wipe dry - in particular when buffer entered the ball-bearing head. Grease the two gliding axes (5) and the nipple (6) at the homogenizing unit daily. Check the holding-screws of the motor and the rack and tighten them if necessary. Replace the aluminium bars holding the bags when they look worn. They are made of „soft“ aluminium to prevent the ball-bearing head from damage.

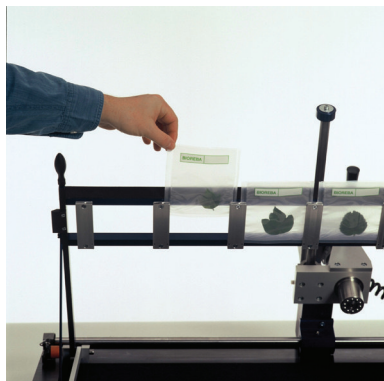
HOMEX 6 contains parts which may rust. Store and operate the HOMEX 6 at a dry place to avoid corrosion. After each use, oil all metal parts with a soft cloth or by using an anti-corrosion protection spray.

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