HOMEX XS

User Guide
 Benutzerhandbuch
 Mode d'emploi

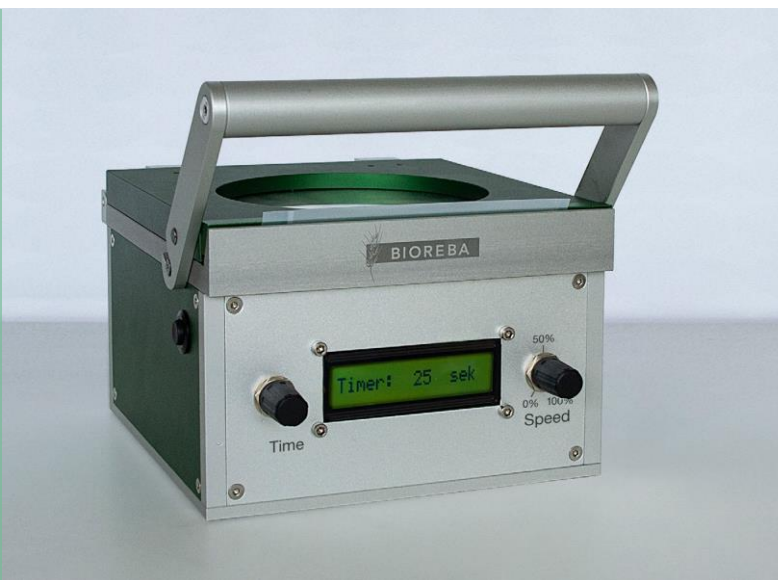
Table of contents / Inhalt / Sommaire

1.	Foreword / Vorwort / Avant-propos	3
1.	Content check / Inhaltsprüfung / Vérification du contenu	3
2.	Safety warnings and guidelines / Sicherheitshinweise und Richtlinien / Consignes et directives de sécurité	3
3.	Maintenance of the instrument / Wartung des Gerätes / Maintenance de l'appareil	5
4.	Introduction / Einführung / Introduction	5
1.	Specifications / Spezifikationen / Spécifications	6
2.	Recommended settings / Empfohlene Einstellungen / Paramètres recommandés	6
5.	Operation guide / Bedienungsanleitung / Guide d'utilisation	7
6.	FAQ	9
7.	Packing list / Stückliste / Conditionnement	10

1. Foreword / Vorwort / Avant-propos

Thank you for purchasing our tissue homogenizer HOMEX XS.

This manual for users describes the instrument's functionalities and contains the operation guidelines. To use the instrument properly, please read this manual carefully before use. Keep it for later reference and troubleshooting.

1.1 Content check

Please check the content according to the packing list (see chapter 7) when you first open the instrument's packaging. If you think there is something wrong with the instrument or something is missing, please contact us immediately.

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Important operation and safety information:

Read the manual before operating the instrument. Read the guidelines below and follow the necessary safety measures as described.

▲ Safe Operating Procedures:

The operation, maintenance and repair of the instrument should comply with the basic usage guidelines and the warnings listed below. Failure to comply will affect the life span of the instrument and will result in the invalidation of the warranty.

▲ Do not:

- use the instrument for anything other than its intended use.
- carry the instrument by the handle.
- use a different power supply than the one supplied (Art. No. 400028).
- open the lid while the instrument is running.
- use the instrument if the cable to the power supply is damaged.

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▲ This product has been designed for indoor and non-extreme temperature conditions use only

▲ The operator should not open or repair the instrument. Opening the instrument will result in a loss of warranty and may cause serious injury or death. If the instrument is not working properly or is defect, only companies approved by the manufacturer should repair it.

▲ The A.C. power's grounding should be sufficiently stable. The 3-pin plug supplied with HOMEX XS power cable is a safety device that should be matched with a suitable grounded socket.

▲ Always place the instrument on a horizontal, stable, and even surface. The surface must be resistant to heat and fluids.

▲ Any liquid in contact with the machine may cause damage and provide a significant hazard.

▲ The instrument should be placed in a dry, cool, and dust-free environment with no exposure to direct sunlight. The instrument should be used only in a well-ventilated environment, with no corrosive gas or strong disturbing magnetic field, away from central heating and other heat sources.

▲ The main switch is on the rear of the device, push "I" to power on the device, and push "O" to power off the device.

▲ Switch off the instrument when you finished work. Pull out the connector plug when there are long periods of inactivity. Cover the instrument with a cloth or paper to protect it from dust.

▲ Pull the connector plug from the back at once in the following case and contact the vendor:

- Liquid flowing into the instrument.
- Drenched or fire burned.
- Abnormal noise or smell.
- Instrument dropped or outer shell damaged.
- The instrument does not function properly.

▲ Keep the connector plug away from heat and damp. If the supply cord is damaged, it must be replaced by the manufacturer, its service agent, or similarly qualified persons, in order to avoid all risks.

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Power off the instrument and pull the connector plug prior to cleaning and clean the HOMEX XS with a damp cloth.

Do not use corrosive cleaning reagents as they may damage the machine. Remove spills of extraction buffer immediately first with a soft cloth wetted with water, then wipe dry, in particular when buffer enters the ball-bearing head.

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4. Introduction / Einführung / Introduction

The HOMEX XS is a rapid, automated system for the homogenization of different plant tissues. The material (e.g. roots, leaves, seeds) is inserted into a BIOREBA extraction bag (Universal 120 x 150 mm or Universal-U-form 120 x 150 mm) which contains an appropriate amount of extraction buffer. The filtered homogenized extract can then be used for downstream applications such as ELISA, DNA/RNA extraction, etc. The HOMEX XS homogenizes any kind of plant tissues without sharp edges (e.g. stones). Some of the samples must be pretreated before extraction.

Seeds need presoaking, while wood slices and thick samples need to be flattened with a rubber hammer. It is also recommended to crush fleshy tissues before extraction.

Up to 100 samples of the size from 0.1 g to 1 g can be processed per hour. The buffer volume used is usually 2 – 5 ml for BIOREBA Universal extraction bags (Art. No. 430100) and should not exceed 3 ml for BIOREBA Universal-U-form extraction bags (Art. No. 440100). Recommended settings are given in table 4.2.

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Note:

- Do not run the machine without an extraction bag inserted as the ball-bearing head (7) might damage the lid.
- The machine will stop automatically when the lid is opened during a homogenization process or if the start button is pressed a second time.

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Figure 1: HOMEX XS front view / Vorderansicht / vue de face

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Figure 3: HOMEX XS rear view / Rückansicht / vue arrière

6. FAQ

■ Instrument won't start homogenization after pressing the start button.

■ Check that all cables have been connected correctly.

■ A sample bag that is too full may prevent the instrument from working. In this case loose slightly the handle until the ball-bearing head starts to rotate. If the instrument runs correctly, tighten again the lid. Scan the QR code (video 1) on this page to watch a short video tutorial.

■ For certain samples (e.g. seeds) it may be necessary to pre-treat them with a rubber mallet. Loosen both screws as shown in Fig. 4 (does not work for all models) and raise/lower the lid according to your samples. Tighten the screws and proceed with homogenization. Scan the QR code (video 2) below for a video tutorial.

■ Homogenate is spilling out of the extraction bag during the homogenization.

■ Please do not exceed the volumes described in table 4.2. When using BIOREBA extraction bags Universal-U-form (Art. No. 440100), a buffer volume of 3 ml and a sample weight of 0.5 g should not be exceeded. For BIOREBA extraction Universal bags (Art. No. 430100), a buffer volume of 5 ml and a sample weight of 1 g should not be exceeded.

■ The sample is not homogenized correctly and non-homogenized particles are moving to the corners.

■ Consider using BIOREBA extraction bags Universal-U-form (Art. No. 440100) instead of BIOREBA bags Universal (Art. No. 430100). Reducing homogenization speed may help preventing particles from moving to the corners.

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EMC-Testcenter AG
Swiss Accreditation Service SAS (STS 0034)
EMC-TESTCENTER AG